


# **VERIcleen®:** Surface Hygiene Indicator



VERIcleen® Test Kit

"In all cases, the results obtained were consistent and easy to interpret and were attainable within one minute of sampling the surface" [Figure 1]

- Ginny Moore and Chris Griffith, "Hygiene checksnthe next generation," International Food Hygiene, vol.13, no.7

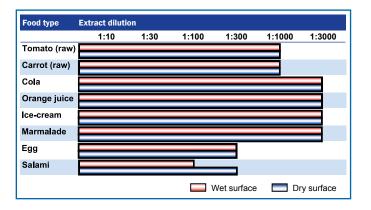

**Figure 1:** The detection of residual food debris from a wet and dry stainless steel surface, using the VERIcleen surface test. Source: Hygiene Checks-the next generation

# **VERICLEEN AT A GLANCE**

- Detects: Organic residue found on unknown surfaces and hand
- Method: Merely wet, wipe, and watch
- >> Economical: Needs no instrument
- ▹ Fast: 5 to 60 seconds
- >> Results: Accurate and repeatable

#### SIMPLIFIED SURFACE MONITORING

VERIcleenTM provides an inexpensive rapid hygiene test with many food/healthcare applications. VERIcleen is simple to use, conveniently packaged and easy to interpret. There is no equipment, no special skills or training needed. VERIcleen is a powerful tool for cleansing programs, hand washing, hygiene compliance, and overall sanitation monitoring.

#### **PROVIDING HYGIENE SAFETY**

Organic residues can be nutrients for bacterial contamination, and they impede disinfectant and sanitizer effectiveness.

VERIcleen validates the quality of cleaning of both product contact surfaces and employee hands by detecting organic material.  $\mathbf{1}$ 

#### **RECOMMENDED FOR:**

Airlines **Auditors Beverage Processors Butcher Shops Caterers Cleaning Companies Commissaries Custom Blending Cruise Ships Dairy Producers Day Care Centers** Delis **Fast Food Chains Food Distributors Food Processors Health Inspectors Hospitals Hotels Nurseries Nursing Homes Restaurants Schools Supermarkets** 

# PROCEDURE



Wet: If surface is dry, wet with 3 or more pumps of Wetting Liquid. Enough liquid must be present on surface to absorb into test when surface is wiped



**Wipe:** Wipe surface with absorbent region of VERIcleen. Check that Wetting Indicator is dark (indicating test strip is sufficiently wet).

NEG/		
	VERIcleen	
POSI	VERIcleen TIVE	
	VERIcleen VERIcleen	

**Watch:** Purple color formation within 1 minute indicates presence of organic residues and need to re-clean.



Supplier Details:

# VERICLEEN: [CODE: VERICLEEN-50K]

- > 50 VERIcleen tests in a black snap cap container
- 1 bottle of Wetting Liquid with atomizer
- Instruction leaflet

### VERICLEEN: [CODE: VERICLEEN-100K]

- >> 100 VERIcleen tests in a black snap cap container
- 2 bottles of Wetting Liquid with atomizer
- Instruction leaflet

